



Waterborne Pathogens Panel

SAMPLE COLLECTION

This panel includes culture for four waterborne pathogens: 1) *Legionella* spp. 2) *Stenotrophomonas maltophilia* 3) *Acinetobacter* spp. 4) *Pseudomonas aeruginosa*.

The sample volume is 500 ml and should be collected as follows:

1. Complete the Chain of Custody form.
2. Use waterproof pen to label bottle with sample location, description, and date.
3. Turn on hot water and immediately fill the bottle (500 ml). Close bottle and invert to mix neutralizer.

SHIPPING

1. Ship the same day samples are collected for receipt in laboratory. The time from sample collection to analysis **should not exceed 30 hours**.
2. Place bottles in box with insulated liner.
3. Place completed Chain of Custody in plastic document holder and pack in box.
4. Ship overnight Monday–Thursday only to:
Special Pathogens Laboratory
1401 Forbes Ave., Suite 401
Pittsburgh, PA 15219

Questions? Contact SPL at 412-281-5335